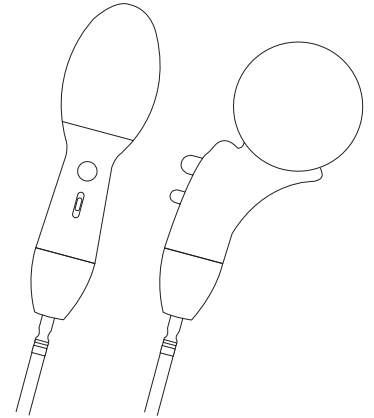
# PROFILLER™ 435/436

## MANUAL PIPETTE CONTROLLERS MANUELLE PIPETTIERHILFEN PIPETEURS MANUELS

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#### SAFETY PRECAUTIONS

- Read operating instructions carefully.
- Observe manufacturer's instructions and guidelines for operation and maintenance.
- Refer to and follow regulations about handling of potentially hazardous reagents.
- Before use, check instrument for good working conditions.
- Do not use harsh chemicals to clean instrument.
- Do not use Profiller™ pipette controllers with liquids incompatible with silicone, PP and ABS, ASA, SAN copolymers.
- Do not allow liquids to enter the instruments.

#### DESCRIPTION

Profiler™ are easy to use manual pipette controllers, which facilitate aspiration and distribution of a variety of reagents. Both instruments are compatible with glass or plastic pipettes with nominal volumes ranging from 1 to 100 mL and perfectly suitable for left or right handed operation.

#### OPERATION

**Pipette fitting:** Squeeze suction bulb. Carefully attach the pipette by inserting it into the holder of the nose cone without using force.

**Aspiration:** hold instrument in a vertical position with the pipette tip down. Place tip of the pipette into the liquid to be aspirated. Press lever control button slowly upward allowing the liquid to fill the pipette until the meniscus level is slightly above the required mark. The suction bulb can be squeezed again if additional aspiration is needed during pipetting. Adjust meniscus position by pressing down the lever control button.

**Dispensing:** place pipette against wall of the receiving vessel and press down lever control button to dispense.

**"To deliver" pipette:** when meniscus has reached a still stand, wait for the required time indicated on the pipette (i.e. EX + 10 s) for dispensing to be completed.

**"Blow out" pipette:** as soon as the meniscus has reached a still stand, press discharge bulb to expel the rest of the liquid contained in the pipette.

**Remove the pipette:** hold the pipette and rotate gently pulling the pipette from the pipette holder.

#### TIGHTNESS

Take care of the instrument to keep it in perfect working condition. To test tightness, attach a pipette and fill it to half of its nominal capacity. Drops falling from the pipette tip would indicate a leak. Unscrew lower assembly and check pipette holder and membrane filter. Reassemble correctly.

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## MAINTENANCE

Do not allow liquid to enter the Profiller™. The PTFE membrane of the protection filter would clog and suction be impaired. To replace the filter, remove the pipette, unscrew nose cone, pull out and replace membrane filter. Choose filter size (0.2 µm or 0.45 µm) according to specific needs. Reassemble the unit. Clean the instrument with a damp cloth. Profiller™ are not autoclavable, except silicone pipette holder.

## WARRANTY

The Profiller™ are guaranteed for a one-year period against any factory or material defects. Damages due to non-respect of manufacturer's instructions are excluded from the warranty. Return the instrument to the dealer from whom it was purchased, after obtaining return authorisation, should regular maintenance not eliminate a detected defect.

**Decontaminate instrument prior to returning it.**

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