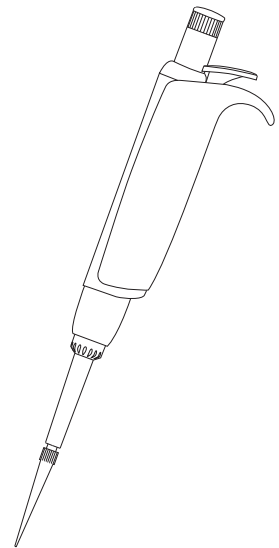
# ACURA<sup>®</sup> manual 826 XS

## EXTRA SHARP MICROPIPETTE MIKROPIPETTE EXTRA SHARP MICROPIPETTE HAUTE PERFORMANCE

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#### SAFETY PRECAUTIONS

- Read operating instructions carefully; keep available for future reference.
- Observe manufacturer's instructions and guidelines for operation and maintenance.
- Refer to and follow regulations about handling of potentially hazardous reagents.
- Before use, check instrument for good working conditions.
- Do not use harsh chemicals (such as acetone) to clean instrument.
- This product should be used only for its intended purpose.
- Mind possible hand-fatigue during serial pipetting and its medical consequences (such as repetitive strain injuries RSI).

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#### DESCRIPTION (Fig. 1)

- Precise and easy pipetting cover volumes ranging from 0.1 µL to 1000 µL.
- Ergonomics and pipette activation provide for an ultra-soft operation.
- Excellent comfort, the short length allows an optimal handling.
- Materials have been carefully selected to ensure durability and optimal resistance to aggressive media.
- Instrument autoclavable at 121°C fully assembled.
- Well positioned, the digital display window (E) allows excellent volume visibility.
- Innovative, integrated and adjustable Justip™ (H) tip ejector for ease of tip ejection.
- Swift-set calibration performed by laboratory personnel (F).

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#### OPERATION (Fig. 1)

##### Holding the pipette

Place finger rest (D) on the phalanx of forefinger. Ease of access of both plunger button (B) and ejector push button (C) allow natural and comfortable thumb movement.

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#### Volume setting

Set desired volume by rotating push button (B) clockwise to decrease the volume and vice versa, until it appears in display window (E). Set volume is indicated by the red index pointer. Precise click-stops of the micrometric mechanism and a free rotating *smartie* colour cap (A) on top of button prevent any volume alteration during pipetting.

**Note:** If the letter E appears beside digits in the display window, the set volume is no longer within working range of the pipette. Overturning the micrometric screw may impair the volume setting mechanism.

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## Tip selection

Disposable Qualitips® - with or without filter - were carefully designed to provide the best pipette and tip combination for all types of applications. Regular tips are autoclavable at 121°C. Filter tips are either autoclavable or supplied in a sterilized version. Use tips supplied by instrument manufacturer, or similar (see tip selection on page 7). Detailed flyer available on request.

Apply new tip as often as deemed necessary in regard with change of sample nature, droplet accumulation after blow out, or foam formation. When using tip racks, place the instrument on tips, press firmly to ensure a good tightness.

**Note:** Enhance precision by rinsing new tip 2-3 times before sampling.

## Tip ejection (Fig. 2)

The patented Justip™ ejection system allows instant height adjustment ( $\pm 2$  mm) of ejector for best fitting of tip used (optimal distance). Set correct position ( $\sim 0.5$  mm space between tip and ejector) by rotating ejector screw ( $< LO - HI >$ ). Click stops prevent any unwanted change while pipetting.

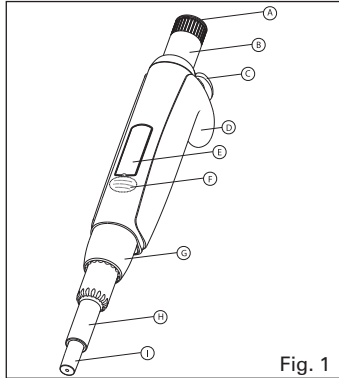
Fig. 1



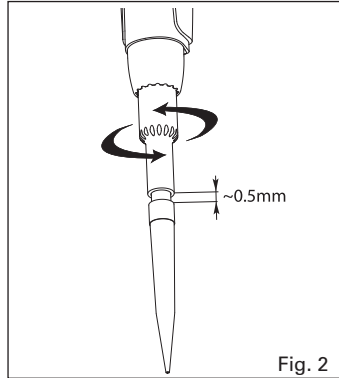
Fig. 2

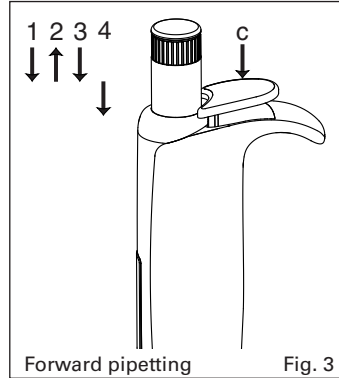Forward pipetting Fig. 3

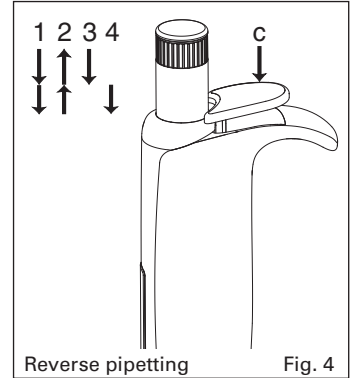Reverse pipetting Fig. 4

## PIPETTING OPERATION

Before pipetting, tightly fit new, clean tips on the nozzle.

### Forward pipetting (Fig. 3)

Exact volume desired is aspirated, then dispensed  
Phase 1 Depress plunger button to intermediate stop ending calibrated stroke.

Phase 2 Immerse tips vertically (2 - 3 mm) into liquid, wait 2 sec. and smoothly release plunger button. Withdraw tips from liquid, without touching container wall.

Phase 3 Place tips against wall of receiving vessel, depress plunger button to first stop and let liquid out.

Phase 4 Depress button to second stop: blow out will expel remaining droplets.

Tip ejection Remove used tips by pressing ejector button (C). Change tips as required.

### Reverse pipetting (Fig. 4)

Liquid is aspirated in excess of the selected volume but only the set volume is dispensed.

Phase 1 Depress plunger button in full.

Phase 2 Immerse tips vertically (2 - 3 mm) into liquid, wait 2 sec. and smoothly release plunger button. Withdraw tips from liquid, without touching container wall.

Phase 3 Place tips against wall of receiving vessel, depress plunger button to first stop only and let liquid out.

Phase 4 Withdraw tips from receiving vessel, sliding it over 10-15 mm along container wall and start operation to phase 2 again, or discard residual liquid by depressing button to second stop.

Tip ejection Remove used tips by pressing ejector button (C). Change tips as required.

## Measuring unknown volumes

The fine volume resolution of the micrometric mechanism allows for precise liquid measurement by aspiration. Remember set volume before starting, dip pipette tip into liquid and turn adjustment button counter-clockwise until unknown quantity is completely aspirated. Measured volume consists in read volume minus initial set volume.

**Example:** volume before aspiration: 10.0  $\mu$ L. Volume read-out after aspiration: 64.7  $\mu$ L. Liquid aspirated: 64.7 - 10.0 = 54.7  $\mu$ L

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## MAINTENANCE

The Acura® manual XS pipettes are in conformity with Council directive IVD 98/79 EEC (in vitro diagnostic medical devices). Consequently, any change on the pipettes (i.e. parts, calibration) is done under responsibility of the user. The models were designed to sustain long-term operation with minimal maintenance and service requirements. However it is recommended to control instrument performance in accordance with laboratory internal procedures (SOP, GLP, etc) or at least once a year. Defect parts must be replaced, or instrument returned to dealer for repair. Order original spare parts from authorised dealers only.

### Tightness control

A tight air chamber is paramount for the instrument to provide accurate and precise results. Leaks can be detected by observing a droplet forming and falling from the extremity of the tips.

Special attention must be paid to the following points in case of leakage:

- Matching of tips and nozzle ?
- Damaged nozzle ?
- Worn-out PTFE sleeve and/or O-ring ?
- Lack of grease ?

### Cleaning

Elements of volumetric module (lower assembly) can be cleaned, or soaked in appropriate decontamination or disinfecting solutions without disassembling.

This operation is particularly important if liquid accidentally entered the pipette. Rinse and dry instrument before reassembling.

An ultra-sonic bath is an efficient way to remove sticking residues, slightly grease O-ring or plunger (see below) before reassembling.

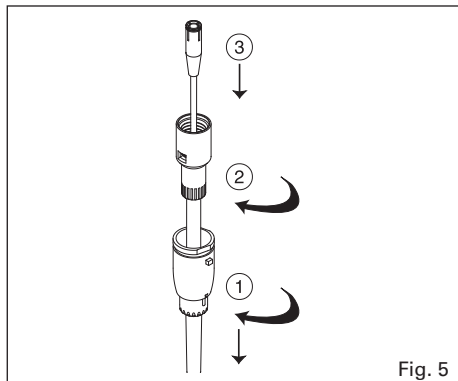
Fig. 5

### Disassembling volumetric module and O-ring change (Fig. 5 and 6)

Removing volumetric module (lower assembly):

- 1) Press ejector button (2), turn part (20) to the left and pull.
- 2) Unscrew barrel (18).
- 3) Depress plunger button in full then unclip (pull) plunger (13).

Barrel assembly (50 - 100 µL):

- To guarantee tightness, minimum friction and spare parts compatibility, the PTFE sleeve can not be changed by itself. Changing the barrel assembly, incl. plunger is necessary.
- Slightly grease plunger (13), sleeve and O-ring and reassemble volumetric module.

Lip seal replacement (200-1000 µL) Fig. 6:

If the instrument is leaking after cleaning (without removing the cylinder head), the following operation is necessary:

- After remove the lower assembly:
  - 1) Lift both clips of the cylinder head (21).
  - 2) Remove the cylinder head (21).
  - 3) Gently remove the lip seal with fingers or with a tip.
- Change the defective parts, clean the plunger (13) and apply thin layer of grease on its surface.
- Slightly grease the lip seal on the top external diameter and between the lips.
- Reposition the lip seal inside the cylinder (18) and clip the cylinder head (21).
- Reassemble the lower assembly.
- When pipetting aggressive solvent, the lip seal (FPM) could get damaged.

**Note:** Tightness parts are not accessible on 2, 10, 10Y and 20 µL models. In case of tightness problem, the whole barrel must be changed. Do not force the plunger into the cylinder.

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## TROUBLE SHOOTING

Observation	Possible cause	Action
Tips not fitting the nozzle	Inadequate position of tip ejector	Adjust position of tip ejector
Inadequate tips	Use original or other compatible tips	
Plunger sticking or dragging	Dirty plunger	Disassemble and clean plunger
Liquid not aspirated	Dry residues block pipette nozzle	Disassemble and clean
Volumetric module not properly assembled	Reassemble according to instructions	
Poor instrument performance	Lack of tightness	Check tip fitting
	Check for damaged nozzle or tightness parts	
Instrument out of calibration	Perform new calibration	
Instrument used with viscous or volatile liquids or at particular temperatures		Perform new calibration with liquid or at current temperature

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### Sterilization

Instrument has been designed for repeated sterilisation in the autoclave at 121°C (20 minutes), fully assembled. Place instrument horizontally in the autoclave, avoiding any direct contact with metal.

Allow instrument to dry and cool down before use. Check tightness and accuracy regularly but at least after 50 autoclaving cycles. Check volumetric module, tighten barrel if loose. Repeated autoclaving may affect material colours and free rotation of *smartie* colour caps. Correct autoclaving and resulting sterility are the responsibility of the user.

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## CALIBRATION

Each Acura® *manual* XS has been factory calibrated and individually controlled. The QC certificate accompanying the instrument includes control data, serial number and operator identification. Calibration is carried out gravimetrically with aqua-bidest. at constant temperature ( $\pm 0.5^\circ\text{C}$ ) comprised between 20 and 25 °C according to ISO 8655.

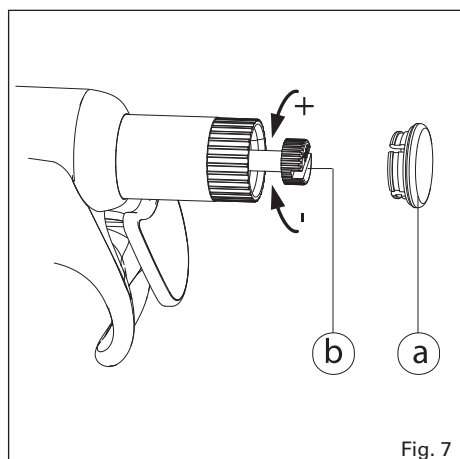
Fig. 7

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### In-lab calibration (Fig. 7 and 8)

The *swift-set* calibration system with integrated key and locking mechanism makes tedious procedures obsolete. If performance results are no longer within recommended values, for instance after QC check or replacing parts, or if changes occur in physical parameters (liquid density, temperature), recalibration is easily performed using an analytical balance. Ten measurements should be performed on small volume after checking seal tightness and perfect working condition of the instrument.

- Remove colour cap (a) on plunger button and pull out calibration screw (b).
- Remove "Calibration seal" (c) and slide calibration cursor (d) on *cal* position (use pointed item such as tip or clip).
- Turn calibration screw until average weighed volume appears in the display window, without holding push button.
- Place calibration cursor on *lock* position. Cover plunger button with *Smartie* cap again.
- Perform few plunger movements and check resulting volume.
- Place a new "Calibration seal" over the calibration cursor.

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## WARRANTY

Your Acura® *manual* XS pipette is guaranteed against any material or manufacturing defects for the period of time specified in its QC certificate. Damages due to non-respect of manufacturer's instructions, safety precautions or autoclaving conditions, as well as material colour alteration are excluded from the warranty. Repair and replacement of parts do not extend warranty time. Should regular maintenance not eliminate a detected defect, return the instrument to the dealer from whom it was purchased after obtaining return authorisation.

**Decontaminate instrument prior to returning it.**

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
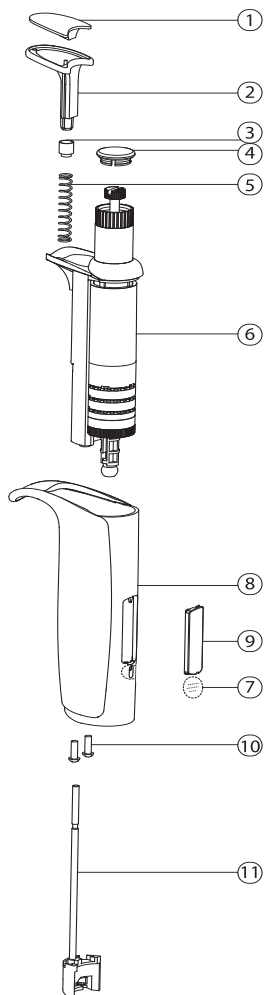
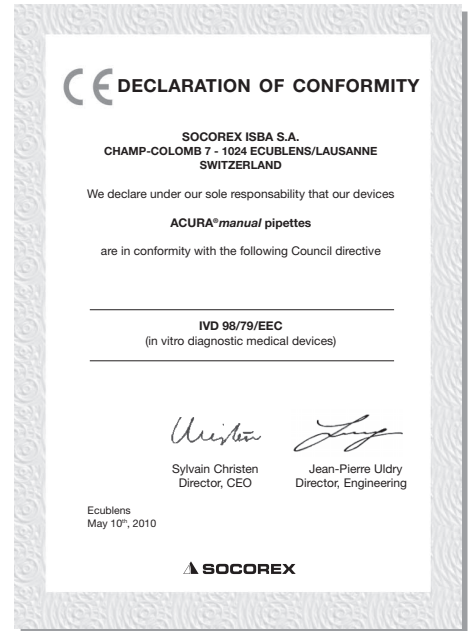
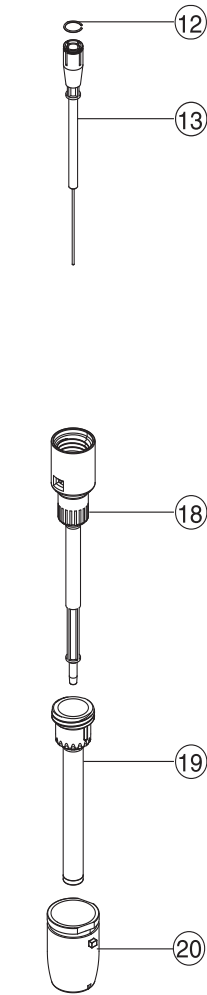
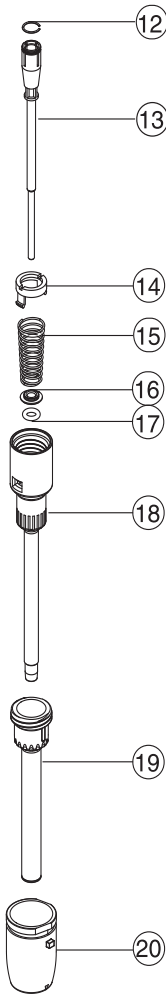
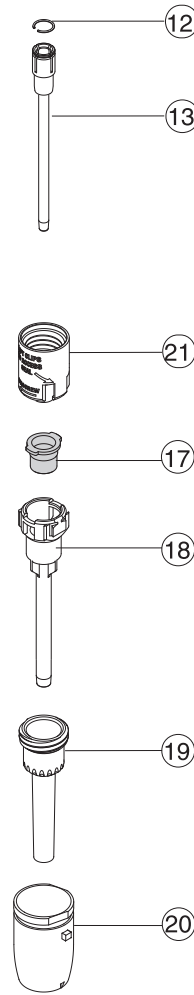
200 µL

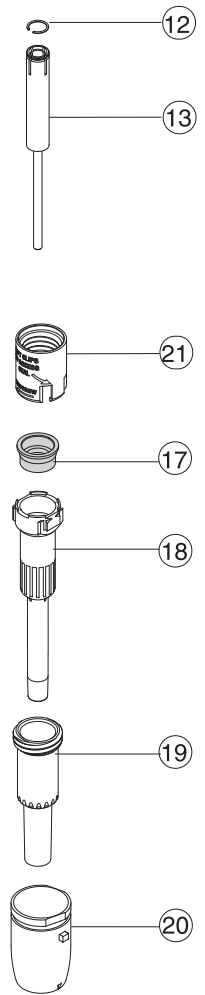1000 µL

**PERFORMANCE AND ORDERING INFORMATION**

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Performance values obtained in forward pipetting mode with bi-dest. water at constant temperature ( $\pm 0.5^\circ\text{C}$ ) between 20 and 25°C in accordance with EN ISO 8655.

**ACCESSORIES**

**ZUBEHÖR**

**ACCESSOIRES**

Description	Packaging	7 positions Model 337	6 positions Model 336
<b>Universal work station – fits single channel pipettes of most brands</b>			
Grey	1 / pack	320.337G	320.336Q
Blue	1 / pack	320.337B	320.336B
Yellow	1 / pack	320.337Y	320.336Y
Green	1 / pack	320.337M	320.336G
Pink / Red	1 / pack	320.337R	320.336R
Orange	1 / pack		320.336O
White	1 / pack		320.336W

Description	Packaging	Socorex Cat. No.
Work station 340 for 3 multi- or single-channel pipettes	1 / pack	320.340
Shelf pipette holder for 2 instruments (transparent blue)	1 / pack	320.332
	4 / pack	320.332.4
Protective calibration seal stickers	25 / bag	1.825.025
	100 / bag	1.825.100

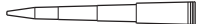
Colour	Cat. No.
<i>Smartie</i> plunger caps, alternative colour selection (6 / pack)	
White	825.710
Grey	825.715
Vanilla	825.716
Lemon	825.712
Yellow	825.717
Orange	825.718
Red	825.714
Rose	825.719
Pink	825.720
Purple	825.721
Ice blue	825.722
Blue	825.713
Mint	825.723
Green	825.711
<i>Smartie</i> mix (14 / pack)	
Assorted colours	1.825.700

**QUALITIPS® PIPETTETIPS**

**QUALITIPS® PIETTENSPIZTEN**

**EMBOUTS QUALITIPS®**

Superior pipette tips cover a range between 0.1  $\mu\text{L}$  and 1000  $\mu\text{L}$ . All models autoclavable at 121°C.

Tips without filter	Description	Vol. ( $\mu\text{L}$ )	Racks	Cat. No.	Bags	Cat. No.
	Nat. ultra-microtip	10	10 x 96/rack	<b>309.0010R</b>	1 x 1000/bag	<b>309.0010B</b>
	Nat. gel load tip	10	4 x 204/rack	<b>302.0010GR</b>		
	Nat. microtip	10	10 x 96/rack	<b>302.0020R</b>	1 x 1000/bag	<b>302.0020B</b>
	Nat. universal tip	200	10 x 96/rack	<b>307.0200R</b>	2 x 500/bag	<b>307.0200B</b>
	Yellow univ. tip	200	10 x 96/rack	<b>327.0200R</b>	2 x 500/bag	<b>327.0200B</b>
	Superior nat. tip	200	10 x 96/rack	<b>308.0200R</b>	1 x 1000/bag	<b>308.0200B</b>
	Nat. gel load tip	200	4 x 204/rack	<b>308.0200GR</b>		
	Nat. tip	200	10 x 96/rack	<b>309.0200R</b>	2 x 500/bag	<b>309.0200B</b>
	Yellow tip	200	10 x 96/rack	<b>329.0200R</b>	2 x 500/bag	<b>329.0200B</b>
	Nat. universal tip	1000	10 x 60/rack	<b>307.1000R</b>	4 x 250/bag	<b>307.1000B</b>
	Blue tip	1000	8 x 60/rack	<b>318.1000R</b>	4 x 250/bag	<b>318.1000B</b>
	Nat. tip	1000	8 x 60/rack	<b>309.1000R</b>	4 x 250/bag	<b>309.1000B</b>
	Blue tip	1000	8 x 60/rack	<b>319.1000R</b>	4 x 250/bag	<b>319.1000B</b>

Efficient protection against aerosols. Tip include high density PE filter for complete, uncontaminated sample retrieval. Certified DNase and RNase free. Sterilized tips are guaranteed pyrogenfree.

Filters tips	Description	Vol. ( $\mu\text{L}$ )	Racks	Cat. No.	Bags	Cat. No.
	Sterilized ultra-microtip	10	10 x 96/rack	<b>309.0010FR</b>		
	Sterilized microtip	10	10 x 100/rack	<b>302.0010FR</b>		
	Autocl. microtip	10			1 x 1000/bag	<b>302.0010AFB</b>
	Sterilized tip	20	10 x 96/rack	<b>308.0020FR</b>	5 x 100/pack	<b>308.0020FS</b>
	Sterilized gel load tip	20	4 x 204/rack	<b>308.0020GFR</b>		
	Sterilized tip	100	10 x 96/rack	<b>308.0100FR</b>		
	Sterilized gel load tip	100	4 x 204/rack	<b>308.0100GFR</b>		
	Sterilized ext. tip	200	8 x 204/rack	<b>308.0200LFR</b>	5 x 100/bag	<b>308.0200LFS</b>
	Sterilized tip	300	10 x 96/rack	<b>308.0300FR</b>		
	Autoclavable tip	300			1 x 1000/bag	<b>308.0300AFB</b>
	Sterilized tip	1000	10 x 100/rack	<b>309.1000FR</b>	5 x 100/pack	<b>309.1000FS</b>
	Autoclavable tip	1000			1 x 1000/bag	<b>309.1000AFB</b>

B = bag, R = rack, F = filter, G = gel, L = long, S = sterilized, A = autoclavable

NOTES

NOTIZEN

NOTES

SOCOREX ISBA S.A. - CHAMP-COLOMB 7 - CH-1024 ECUBLENS/LAUSANNE - SWITZERLAND  
socorex @ socorex.com - www.socorex.com - TEL. +41 21 651 6000 - FAX +41 21 651 6001

SAS Accredited  
Service Laboratory  
ISO/IEC 17025

**G.M. SYSTEM**  
**ISO 9001/13485**  
**CERTIFIED**