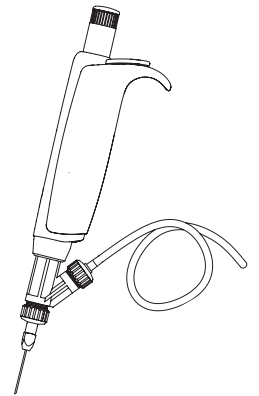
# ACURA® *manual* 865

## MICRODISPENSER PIPETTE MIKRODISPENSIER PIPETTE PIPETTE MICRODOSEUR

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#### SAFETY PRECAUTIONS

- Read operating instructions carefully, keep available for future reference
- Observe manufacturer's instructions and guidelines for operation and maintenance
- Refer to and follow regulations about handling of potentially hazardous reagents
- Before use, check instrument for good working conditions
- Prime and rinse carefully when left unused for a while or before disassembling
- Do not use harsh chemicals (such as acetone) to clean instrument
- This product should be used only for its intended purpose
- Mind possible hand-fatigue during serial pipetting and its medical consequences (such as repetitive strain injuries RSI)

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#### DESCRIPTION

The Acura® *manual* 865 is a hand-held microdispenser pipette intended for repetitive, precise distribution of volumes ranging from 5 to 1000 µL. Based on the Acura® *manual* 825 model, it includes a three-way valve system for automatic self-refilling. After priming, the instrument operates without air interface in the barrel chamber.

Major features include large display window for permanent volume visibility, superior ergonomics and ultra-soft operation. Calibration easily performed by laboratory personal thanks to *swift-set* system.

Carefully selected materials ensure durability and optimal resistance to aggressive media. The unit is autoclavable at 121°C fully assembled.

Each microdispenser pipette is supplied with a 90 cm silicone tubing, blunt end cannula (0.8 × 40 mm), female Luer inlet valve cap, QC certificate and operating instructions.

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## OPERATION

### Holding the microdispenser pipette

Place finger rest on the phalanx of forefinger. Ease of access of plunger button allows natural and comfortable thumb movement. The valve system revolves for optimal feed source positioning.

### Volume setting

Set desired volume by rotating push button clockwise to decrease the volume and vice versa, until it appears in display window. Selected volume is indicated by the top of the pointer on window. Precise click-stops of the micrometric mechanism and a free rotating *smartie* colour cap on top of button prevent any volume alteration during dispensing.

**Warning:** When setting a smaller volume on a primed dispenser pipette, liquid in the barrel will be expelled through the outlet valve.

**Note:** If the letter E appears beside digits in the display window, the set volume is no longer within working range of the pipette. Overturning the micrometric screw may impair the volume setting mechanism.

### Cannula selection (Fig. 1 and 2)

One stainless steel blunt end cannula (0.8 × 40 mm) is included (1). Additional sizes are available on request. Select cannula size in relation with the distributed liquids. A too small internal diameter make plunger activation harder and modify instrument performance.

To fit the optional PTFE cannula (2), remove outlet valve cap (No. 22) and tightly screw PTFE cannula in place.

Use the optional manifolds (4 or 8 ×) for multichannel distribution. Use of Luer Lock outlet valve cap is recommended for better fitting of manifold.

**Note:** divide set volume by number of channels to obtain volume distributed per channel.

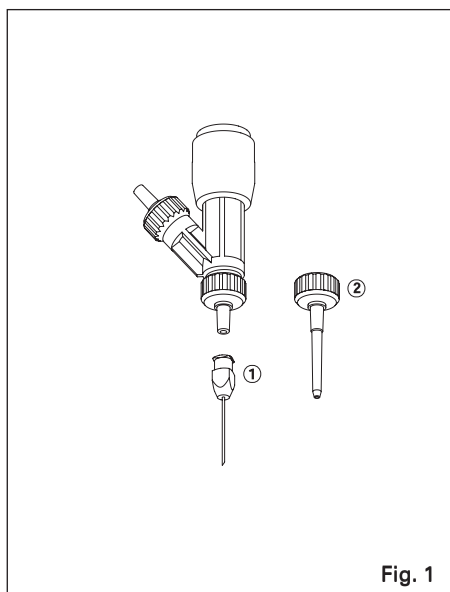
Fig. 1

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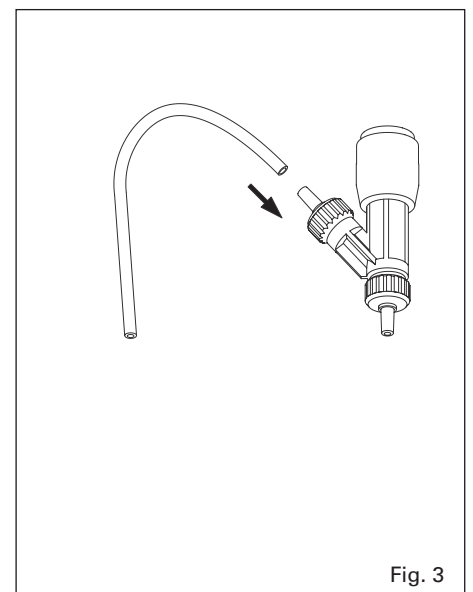
Fig. 3

### Feed source selection (Fig. 3)

Connect 90 cm silicone tubing supplied to male Luer intake valve cap (No. 22).

Alternative sources:

Fill in and connect 30 mL PP or PTFE reservoirs directly to male Luer intake valve cap, then prime microdispenser (Fig. 4).

When using all-glass or plastic disposable syringes (Fig. 5), change intake valve cap for female luer cap (cat. No. 861.720). Fill in syringe then connect to intake valve cap. Prime before dispensing.

For vaccine-type vial (Fig. 6), puncture rubber stopper with feed needle of optional vial holder (1) and slide bottle lock into place (2). Fix to male Luer intake valve cap (3). Prime before dispensing.

### Priming

Activate several times plunger fully set on nominal volume to replace air with liquid. Primed valuable liquid should be collected in a clean vessel for recycling. Insufficient priming may alter instrument performance.

### Dispensing

Press push-button to distribute liquid. Constant aspirating / dosing speed provides for better result consistency. Never apply force. Adapt dosing speed to avoid splashing.

### End of dosing

When dosing is completed, remove feed source. Empty barrel chamber and valves completely by pressing push-button. Flush valves with appropriate detergent and thoroughly rinse with distilled water.

## MAINTENANCE

The microdispenser pipette was designed to sustain long-term operation with minimal maintenance and service requirements. However it is recommended to control instrument performance in accordance with laboratory internal procedures (SOP, GLP, etc.) or at least once a year. Defect parts must be replaced, or instrument returned to dealer for repair. Order original spare parts from authorised dealers only.

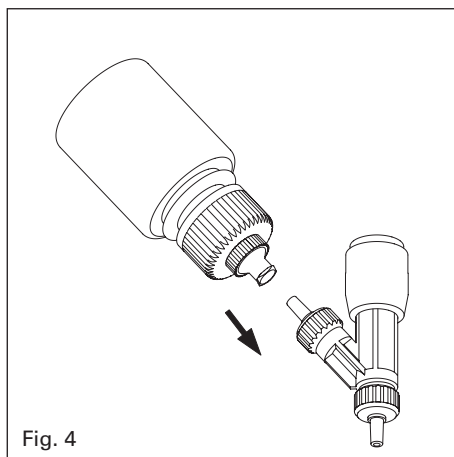
Fig. 4

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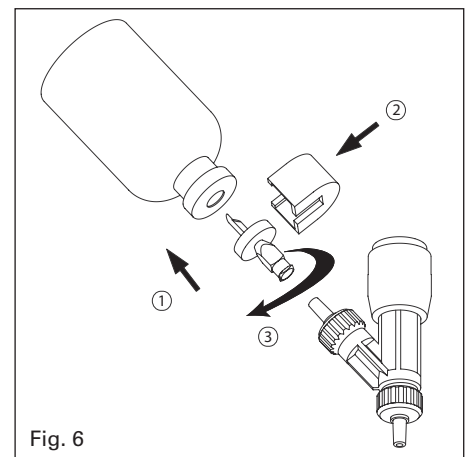
Fig. 6

A clogged aspiration valve can easily be fixed up by detaching the ball from the valve seat. Therefore, introduce a small rod into the intake valve opening and gently push until the ball moves.

**Note:** using force may damage valve spring.

### Tightness control

Perfect tightness is paramount for accurate and precise results. Intake and outlet valves should be tightly screwed to ensure optimal dosing. Particles in suspension may impair valve function. If necessary, liquid should be filtrated prior to distribution.

### Cleaning

Plunger and valve assembly can be cleaned with alcohol, or soaked in appropriate detergent or disinfecting solutions.

If ultra-sonic bath is used to remove sticking residues in barrel, slightly grease O-ring before reassembling.

After long activation, some liquid may be found in the back of the plunger. When dosing is completed, remove valve unit and clean plunger. Check proper tightness after reassembling.

### Disassembling – reassembling (Fig. 7-9)

Unscrew valve body (No. 19). Depress plunger button in full then unclip (pull) plunger (No. 14).

Remove valve caps (No. 22) to access to the valves. Remove barrel sleeve (only on 50 and 200 µL models) and clean barrel chamber.

Plunger and PTFE sleeve do not need to be greased. Grease O-ring slightly if sleeve was disassembled.

Reassemble plunger and parts of valve body according to figure 9. Aspiration (No. 21) and dosing (No. 20) valves have different shapes to prevent mispositionning.

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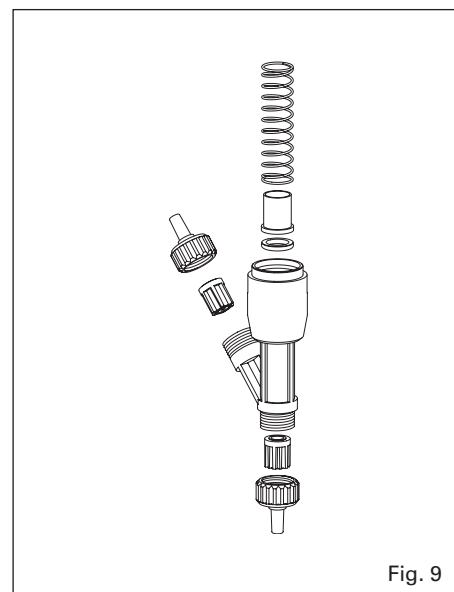
Fig. 9

## Materials

Parts in contact with liquid are chemically inert:

- Valve body: PVDF
- Valve seat: sapphire glass
- Valve ball: synthetic ruby
- Valve spring: platinum-iridium
- Valve cap: PVDF
- Barrel assembly: PVDF
- Plunger: inox DIN 316L (50 µL) or PVDF
- Plunger sleeve: PTFE
- Plunger O-ring: FPM

**Warning:** Damages due to incompatible reagents are not warranty covered. In case of doubt, check for chemical compatibility (see below) or refer to manufacturer before use.

## Sterilization

Instrument has been designed for repeated sterilisation in the autoclave at 121°C (20 minutes), fully assembled. Place instrument horizontally in the autoclave, avoiding any direct contact with metal. Allow instrument to dry and cool down before use. Check tightness and accuracy regularly but at least after 50 autoclaving cycles. Check valve assembly, tighten lower assembly and valve caps if loose. Repeated autoclaving may affect material colours and free rotation of *smartie* colour caps. Correct autoclaving and resulting sterility are the responsibility of the user.

## CALIBRATION

Each microdispenser pipette has been factory calibrated and individually controlled. The QC certificate accompanying the instrument includes control data, serial number and operator identification. Calibration is carried out gravimetrically with aqua-bidest. at constant temperature ( $\pm 0.5^\circ\text{C}$ ) comprised between 20 and 25 °C according to ISO 8655, parts 5 and 6.

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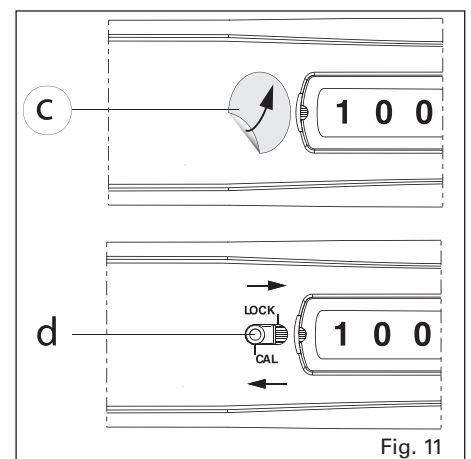
Fig. 11

### In-lab calibration (Fig 10-11)

The *swift-set* calibration system with integrated key and locking mechanism makes tedious procedures obsolete. If performance results are no longer within recommended values, for instance after QC check or replacing parts, or if changes occur in physical parameters (liquid density, temperature), recalibration is easily performed using an analytical balance. Ten measurements should be performed on small volume after checking seal tightness and perfect working condition of the instrument.

- Remove colour cap (a) on plunger button and pull out calibration screw (b).
- Remove "Calibration seal" (c) and slide calibration cursor (d) on *cal* position (use pointed item such as tip or clip).
- Turn calibration screw until average weighed volume appears in the display window, without holding push button.
- Place calibration cursor on *lock* position. Cover plunger button with *Smartie* cap again.
- Perform few plunger movements and check resulting volume.
- Place a new "Calibration seal" over the calibration cursor.

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### WARRANTY

Your Acura® *manual* microdispenser pipette is guaranteed against any material or manufacturing defects for the period of time specified in its QC certificate. Damages due to non-respect of manufacturer's instructions, safety precautions or autoclaving conditions, as well as material colour alteration are excluded from the warranty. Repair and replacement of parts do not extend warranty time.

Should regular maintenance not eliminate a detected defect, return the instrument to the dealer from whom it was purchased after obtaining return authorisation.

**Decontaminate instrument prior to returning it.**

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### PERFORMANCE AND ORDERING INFORMATION

| Volume range (µL) | Division (µL) | Inaccuracy* (Err.Rel%) |           | Imprecision* (CV%) |           | Cat. No. |
|-------------------|---------------|------------------------|-----------|--------------------|-----------|----------|
|                   |               | Min. vol.              | Max. vol. | Min. vol.          | Max. vol. |          |
| 5 - 50            | 0.1           | < ± 5.0%               | < ± 1.5%  | < 2.0%             | < 0.4%    | 865.0050 |
| 20 - 200          | 0.2           | < ± 2.5%               | < ± 1.0%  | < 1.5%             | < 0.3%    | 865.0200 |
| 100 - 1000        | 1             | < ± 1.5%               | < ± 0.6%  | < 0.6%             | < 0.2%    | 865.1000 |

\*Performance values obtained by a regular and smooth plunger movement, with *Acqua bidest.* at constant temperature (± 0.5°C) comprised between 20 and 25°C

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\*\*Cinematic viscosity indicated in centistokes (cst). Test conditions: bottle feeding, aspiration time < 3 seconds, no cannula. Dynamic viscosity in centipoises (cps) is obtained by multiplying the cinematic viscosity by the density of the liquid (kg/m³).

### ACCESSORIES

### ZUBEHÖR

### ACCESSORIES

| Description   | Packaging | Cat. No.  |
|---|-----------|-----------|
| Reservoir with Luer connector, PP, 30 mL                  | 1 / pack  | 1.861.631 |
| Reservoir with Luer connector, PTFE, 30 mL                | 1 / pack  | 1.861.21T |
| Vial holder with feed needle and lock                     | 1 / pack  | 1.861.925 |
| Inlet valve cap, female Luer (syringe feed)               | 1 / pack  | 1.861.720 |
| Outlet valve cap, Luer lock                               | 1 / pack  | 1.861.569 |
| SS delivery cannula, 0.8 × 40 mm                          | 12 / pack | 370.0840  |
| PTFE delivery cannula                                     | 1 / pack  | 1.861.E32 |
| Work station 340 for 3 pipettes                           | 1 / pack  | 320.340   |
| Shelf pipette holder for 2 instruments (transparent blue) | 1 / pack  | 320.332   |
|   | 4 / pack  | 320.332.4 |

| Description  | Packaging | Cat. No.  |
|--|-----------|-----------|
| <i>Smartie</i> mix, assorted colours   | 14 / pack | 1.825.700 |
| <i>Smartie</i> plunger cap, blue   | 6 / pack  | 1.825.713 |
| <i>Smartie</i> plunger cap, yellow   | 6 / pack  | 1.825.717 |
| <i>Smartie</i> plunger cap, orange   | 6 / pack  | 1.825.718 |
| See Acura® <i>manual</i> pipette line for additional <i>smarties</i> colours |           |           |
| Silicone feed tube, 90 cm  | 1 / pack  | 861.719   |
| 4-channel manifold, Luer lock  | 1 / pack  | 1.170.054 |
| 8-channel manifold, Luer lock  | 1 / pack  | 1.170.058 |
| Protective calibration seal stickers   | 25 / bag  | 1.825.025 |
|  | 100 / bag | 1.825.100 |

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