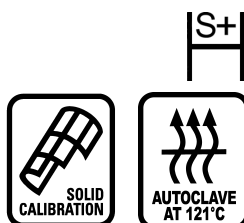
CALIBREX™ 521

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SAFETY PRECAUTIONS

- Read operating instructions carefully, observe manufacturer's general recommendations
- Refer and follow regulations about handling of potentially hazardous reagents
- Before each dosing operation, control perfect working, tightness condition and stability on bottle
- Prime and rinse dispenser carefully when left unused for a while or before disassembling
- Delivery jet should never point towards a person while dosing
- Blocked parts should not be separated by applying force on plunger/barrel assembly or on sleeve
- Remove tip cap from delivery jet before dosing
- Maximal temperature of use < 50°C
- Cover delivery jet with cap and hold reagent reservoir firmly when carrying dispenser
- This instrument should be used only for its intended purpose

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DESCRIPTION

Built with carefully selected materials, the adjustable volume Calibrex dispenser 521 is autoclavable fully assembled. It allows a safe repeated dosing of a large number of reagents with outstanding accuracy and precision. Three models are available to cover a range from 2 to 100 mL (2 – 20 mL, 10 – 50 mL, 10 – 100 mL). Three other models of Calibrex 520 cover smaller volume ranging from 0.25 to 10 mL.

The instrument fits directly most laboratory bottles with standard Ø 32 mm screw neck. Threaded adapters for Ø 38, 40, 45 mm are supplied. Additional diameters and adapters for tapered necks are available upon request.

The Solid Calibration System®, consisting in one cam with pre-calibrated steps provides for digital volume entry and numerical display. An additional fine tuning ring allows quick volume setting. The dispenser Calibrex 521 is easy to operate, maintain and calibrate.

The Calibrex 521 fixed volume models have the same characteristics, but volume is locked on a selected position.

OPERATION

Calibrex 521, 20 mL is supplied fully assembled. On 50 and 100 mL models, first remove shipping plug and attach delivery jet assembly No 13. Introduce without applying force feed tube No 9 in the aspiration valve No 6. Cut feed tube to the desired length if necessary.

Select an adapter appropriate to the bottle and install the instrument. Use a stand to hold the unit if original reagent bottle does not provide for sufficient base stability (Fig 1) see accessories on page 4.

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DOSING INSTRUCTION

Volume setting

The selected volume is shown in the windows of the guide bar No 7 (fig 2). Rotate selector No 2 until desired main volume appears. It automatically clicks-in on the right position. While holding selector, rotate vernier No 16 until desired intermediary volume appears in the windows. Correct position is achieved when both numbers are on one line.

Priming

To prime the unit, activate the plunger fully at least once and follow with several short strokes up to lower stop to eliminate air bubbles in the delivery jet. Liquid should be collected in a clean vessel and may be recycled.

Dispensing

Remove tip cap No 14 before dosing. Raising the plunger fills the dispenser barrel. Depress to dispense liquid, a drop may appear. Best results are obtained by moving the plunger gently and regularly between the upper and the lower stop. Liquid density or viscosity may change dosing speed and instrument performance. Never apply force when working.

End of dosing

After use, place tip cap No 14 on delivery jet to prevent fluid from leaking back, and valves to dry out. Rinse instrument carefully before storing if liquid likely to solidify.

MAINTENANCE

Regular maintenance of the Calibrex dispenser 521 will substantially contribute to its optimal working and long life span.

Opening

Make sure the dispenser barrel contains no liquid by depressing plunger down to lower stop. Remove dispenser from the bottle and prime if necessary. Remove feed tube No 9. Unscrew stop screw No 12 and pull cap No 1 to take out plunger from barrel. Unscrew ring No 19 and pull dispenser body No 5. This allows access to barrel No 4. Remove delivery jet No 13 to access to valve No 10. After reassembling, make sure the dispenser works properly. Damaged parts should be exchanged, using original spare parts only.

Cleaning

Clean fluid path of the dispenser regularly, i.e. whenever the bottle has been emptied or before dosing another liquid.

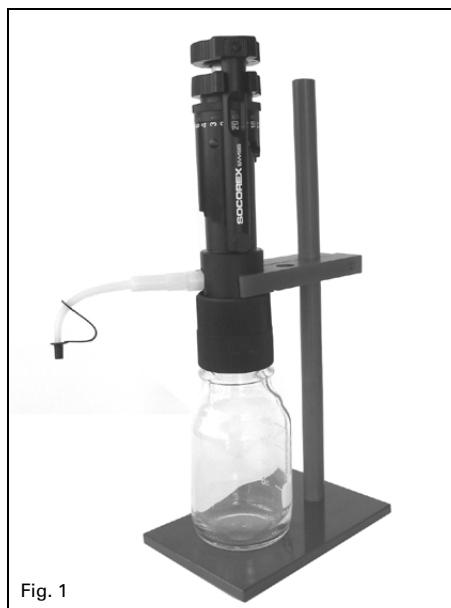
Fig. 1

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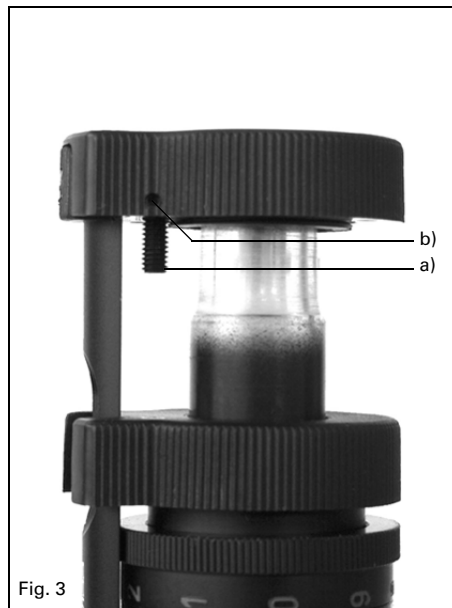
Fig. 3

When plunger is dragging or sticking disassemble the instrument. Soak dirty parts in a mild laboratory detergent. Rinse carefully then check correct function after assembling. To separate a frozen plunger, never use force, the glass barrel could break. Leave frozen plunger assembly soaking in an appropriate detergent overnight. Check regularly both aspiration valve No 6 and dispensing valve No 10. Ultra-sonic bath may help to remove residue on dirty valves.

Rassembling

Damaged parts should be exchanged, using only original spare parts. After reassembling, make sure all connections are tight and check correct function of the instrument.

Calibration

The instrument is factory calibrated and does not require any particular calibration. Should the performance of the instrument during routine control no longer correspond to the manufacturer's specifications or if conditions of use (liquid density, temperature, etc.) require recalibration, this can be easily performed gravimetrically using an analytical balance.

A control and calibration of the dispenser is also recommended when changing following parts: Plunger cap No 1, selector No 2, plunger and barrel assembly No 4, guide bar No 7 or vernier ring No 16.

Calibration screw a), located under plunger cap No 1, corresponds to the lower stop of plunger travel (fig 3). Shorten the lower stop position (calibration screw deeper in cap No 1) will increase the volume and vice versa. Unscrew calibration fixing screw b) before performing a calibration and check resulting volume (smooth dispensing at constant speed) after each adjustment. Tighten fixing screw again after calibration.

Sterilisation

Instrument has been designed for repeated sterilisation in the autoclave at 121°C (20 minutes), fully assembled. Do not sterilize an instrument which has not been primed and rinsed. Place instrument horizontally in the autoclave, avoiding any direct contact with metal. Allow instrument to dry and cool down to room temperature before use. Make sure plunger and all connections are securely tight. Repeated autoclaving may affect material colours. Correct autoclaving and resulting sterility are the responsibility of the user. Calibrex 521.020 may be autoclaved when fitted with the PTFE extension tubing (cat. No 1.524).

Materials

All parts in contact with the liquid are chemically inert:

- Valve body: ETFE / ceramic
- Valve balls: synthetic ruby or ceramic
- Valve seat: glass
- Valve spring: platinum-iridium
- Barrel: neutral borosilicate glass
- Plunger: glass with sturdy PFA coating
- Body: ETFE
- Feed tube and delivery jet: PTFE/ETFE

Warning: Vapours of concentrated strong acids may damage your instrument after some time and reduce its life span. It is recommended not to leave the dispenser permanently on a bottle containing concentrated strong acids and to rinse it after use. To prevent long term loss of product by evaporation, the same recommendation is valid when using volatile solvents.

Damages due to incompatible reagents are not warranty covered. In case of doubt, check for chemical compatibility or refer to manufacturer before use.

WARRANTY

The 521 Calibrex dispenser is guaranteed for a period of one year against any factory or material defect. Glass breakage or cases due to non respect of manufacturer's and safety instructions are excluded from the warranty. Repairs and changing parts do not extend the warranty period.

Should regular maintenance not eliminate a detected defect, return the instrument to the dealer from whom it was purchased, after obtaining return authorization.

Decontaminate instrument prior to returning it.

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PERFORMANCE DATA AND ORDERING INFORMATION

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
*Values refer to the maximum volume, obtained by a regular and smooth plunger movement, with Acqua bidest. at constant temperature ($\pm 0.5^\circ\text{C}$) comprised between 20 and 25°C, in accordance with ISO 8655. ** When ordering, specify desired fixed volume within instrument range.

Vol. (mL)	Max. cinematic viscosity (cst)***	Example of liquids	Test conditions for all 521 models
20	100	Olive oil	Dispensing time < 10 s Dispensing strength < 3 kg Pressure in barrel < 3 bars
50	300	Silicone oil	
100	150	Silicone oil	

***Cinematic viscosity is indicated in centistokes (cst). The dynamic viscosity in centipoises (cps) is obtained by multiplying the cinematic viscosity by the density of the liquid (kg/m^3).

Cat. No	Description	Cat. No	Description	Cat. No	Description
GLP.xx	Threaded adapter, PP, Ø 22, 25, 28, 30, 34, 36, 38, 40, 45 mm	314.xxxx	Amber glass reservoir (Ø 32/45 mm) 0100, 0250, 0500, 1000, 2500 mL	314.0500P	Pyrex glass reservoir (Ø 32 mm) 500 mL with side neck
GLT.xx	Threaded adapter, PTFE, Ø 36, 38, 40, 45 mm	314.xxxxPE	Amber glass reservoir with PE coating (Ø 32/45 mm) 0500, 1000, 2500 mL	315.xxxx	PE reservoir (Ø 25/45 mm) 250, 0500, 1000, 2500 mL
NSP.xx	Tapered adapter, PP, 18.8, 24, 29.2, 45 mm	320.SB050	Stand for one dispenser (up to 50 mL model)	521.621	Spare feed tube 300 mm, Ø 7,5 mm, PTFE, 20 mL model
NST.xx	Tapered adapter, PTFE, 18.8, 24, 29.2 mm	320.BC050	Stand for remote liquid intake (20 and 50 mL models)	521.622	Spare feed tube 300 mm, Ø 10 mm, PTFE, 50 mL model
1.524	60 cm delivery extension tubing, PTFE, for 20 mL model	320.BC100	Stand for remote liquid intake (100 mL model)	521.623	Spare feed tube 375 mm, Ø 10 mm, PTFE, 100 mL model

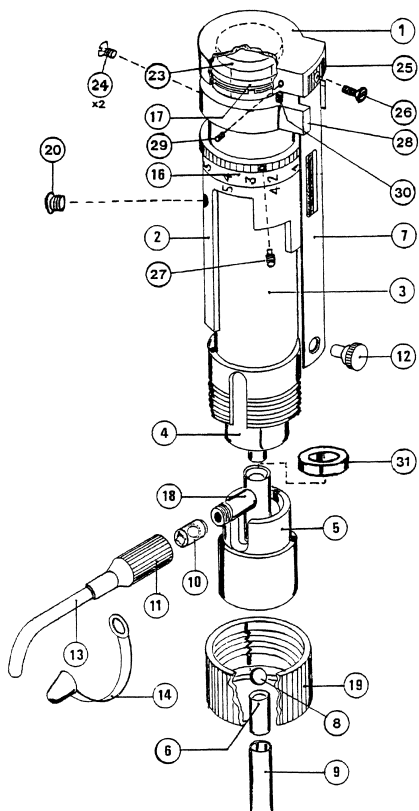SPARE PARTS

- Plunger cap
- Selector assembly
- Sleeve
- Glass barrel
- Body
- Intake valve
- Guide bar
- Intake valve ball
- Feed tube
- Outlet valve assembly
- Delivery jet connector
- Stop screw
- Delivery jet
- Tip cap
- Vernier ring
- Nut
- T-union
- Ring
- Clik stop assembly
- Guide ring fixing screw
- Guide bar fixing plate
- Guide bar fixing screw
- Vernier finger
- Guide ring
- Calibration fixing screw
- Calibration screw
- Distance ring

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