STEPPER™ 411 / ECOSTEP™ SYRINGES

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SAFETY PRECAUTIONS

- Read operating instructions carefully; keep available for future reference
- Observe manufacturer's instructions and guidelines for operation and maintenance
- Refer to and follow regulations about handling of potentially hazardous reagents
- Before use, check instrument for good working conditions
- Use only original Ecostep™ syringes
- Use water or alcohol to clean instrument (do not use solvents such as acetone)
- Mind possible hand fatigue and its medical consequences
- This product should be used only for its intended purpose

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DESCRIPTION

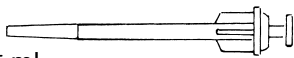
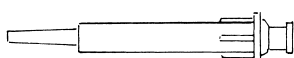
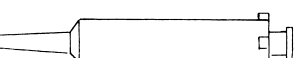
The Stepper™ 411 is a unique trigger action repeater pipette for fast liquid dispensing without thumb fatigue. Careful design provides for comfort and trouble free dispensing of up to 73 samples without refilling. Sample volume and number of shots are clearly displayed on setting knob. Only three sizes of single use Ecostep™ syringes cover from 10 µL to 5 mL. They are the only part in direct contact with the liquid which greatly reduces risks of carry over or cross contaminations. The Ecostep™ syringes are available bulk or as high purity grade, single wrapped sterilized *bioproof*™ certified version. An automatic locking mechanism prevents dispensing errors when the syringe is almost empty.

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Three setting knobs of different colours – one for each syringe size – are supplied with the Stepper. The box includes samples of the Ecostep™ syringes.

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OPERATION

Syringe selection

Non sterile Ecostep™ supplied in bulk are perfectly adapted for all routine applications.

When sterility or high purity grade are required, for microbiology or molecular biology applications, choose the single wrapped sterilized Ecostep *bioproof*™. Each production batch is tested to be sterile and free of detectable human DNA, DNase, RNase and pyrogens (endotoxins). Detection levels indicated in batch certificate. See ordering information p 4.

Made of PE / PP, the Ecostep™ syringes show overall good chemical resistance. Use QR code below for instant access to chemical resistance chart. Damages due to incompatible reagents are not covered by the warranty.

Inserting the syringe and setting button

- Select button (3) and colour matching syringe (5) according to the volume to be delivered.
- Insert button (3) on shaft as shown in picture (B), making sure guide (8) enters slot (9) in button. To remove button, just pull out strongly.
- Slide rod (2) down until it stops.
- Insert syringe (5) into recess (4) until it stops. Red syringe needs adapter (7), supplied with the instrument, to be fitted first (Fig. 1).

Setting the volume (Fig. 2)

- Unlock setting button (3) by pressing it against casing and rotate until selected volume on scale is even with setting mark (6) on casing.
- Release button (3) and allow to snap back neatly into locking position. Together with the delivered volume, the scale shows the number of shots available when syringe has been filled completely.

The syringe graduation helps evaluate the number of shots available when the syringe is partially filled.

DOSING INSTRUCTIONS

Filling the syringe (Fig. 3)

- Immerse nozzle of syringe (5) into liquid.
- Fill by **slowly** pulling rod (2) upwards until it stops or, if syringe is to be partially filled, up to the corresponding graduation mark. Small air bubbles on the plunger will not affect precision.
- Press trigger (1) once until it stops to position stepper mechanism on departure line. **Discard first delivery.** When working with expensive reagents, it may be advisable to switch setting button (3) to smallest volume and to dial the actual sample volume only after the first delivery has been discarded.

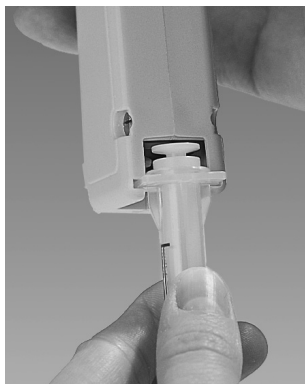
Fig. 1

Fig. 2

Fig. 3

Fig. 4

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Dispensing (Fig. 4)

A yellow pipette tip attached to the yellow and blue syringes will facilitate the access into narrow tubes and increase the precision of the delivery.

- Place syringe against wall of receiving vessel and depress trigger (1) until it stops; release completely when delivery has been completed.
- To prevent dispensing errors, an automatic device locks the trigger whenever the liquid remaining in the syringe at the end of a run is less than the set volume. The residue can be cleared by pushing rod (2) down.
- Dispensing speed should not exceed 1-2 shots/sec.

Removing the Ecostep™ syringe

- Push rod (2) down until it stops.
- Slide syringe out of recess (4).
- Red syringe adapter can be reused.

MAINTENANCE

The Stepper™ 411 does not need any maintenance. Do not open the instrument casing. In case of technical problem, send it to your local dealer for repair after obtaining return authorization. **Decontaminate instrument prior to returning it.**

Calibration

Each Stepper™ is supplied with an individual QC certificate. There is no need and no possibility to calibrate the instrument. Check performance using new, clean Ecostep™ syringes under stabilized temperature condition (20-25°C ± 0.5°C) for instrument, syringe and distilled water. If performance is incorrect, return to dealer for servicing.

Sterilization

The Stepper™ and Ecostep™ syringes are not autoclavable. For sterile applications, order sterilized, single wrapped Ecostep *bioproof*™ syringes.

WARRANTY

The Stepper™ is guaranteed for a period of two years against any factory or material defect. Cases due to non-respect of safety precautions and manufacturer's instructions as well as colour alteration are excluded from the warranty. Repair and replacement of parts do not extend the warranty time.

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Chemical compatibility

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PERFORMANCE DATA AND ORDERING INFORMATION

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Volume	Syringe colour	Syringe capacity	Inaccuracy (E%)	Imprecision (CV%)	Cat. No.
10 – 100 µL	Yellow	0,75 mL	≤± 1.5% at 20 µL ≤± 1.0% at 100 µL	< 1.5% < 0.8%	411.5000
50 – 500 µL	Blue	3,75 mL	≤± 1.0% at 100 µL ≤± 0.5% at 500 µL	< 1.0% < 0.5%	
500 – 5000 µL	Red	27,5 mL	≤± 0.8% at 1000 µL ≤± 0.5% at 5000 µL	< 1.2% < 0.4%	

Performance test in accordance with ISO 8655 method using bi-distilled water at constant temperature (+/- 0.5°C) between 20 and 25°C. Stepper 411 include three colour coded selector buttons, adapter, Ecostep™ syringe samples.

Ecotstep Syringe colour	Version	Packaging	Cat. No.
Yellow	Non-sterile	Bulk, 100/box	316.010
	Sterile <i>bioproof</i> ™ single wrapped	100 x 1/box	316.010.9
Blue	Non-sterile	Bulk, 100/box	316.050
	Sterile <i>bioproof</i> ™ single wrapped	100 x 1/box	316.050.9
Red	Non-sterile	Bulk, 100/box	316.500
	Sterile <i>bioproof</i> ™ single wrapped	100 x 1/box	316.500.9

Ecotstep *bioproof*™ syringes are certified free of detectable human DNA, DNase, RNase and pyrogen.

ACCESSORIES

ZUBEHÖR

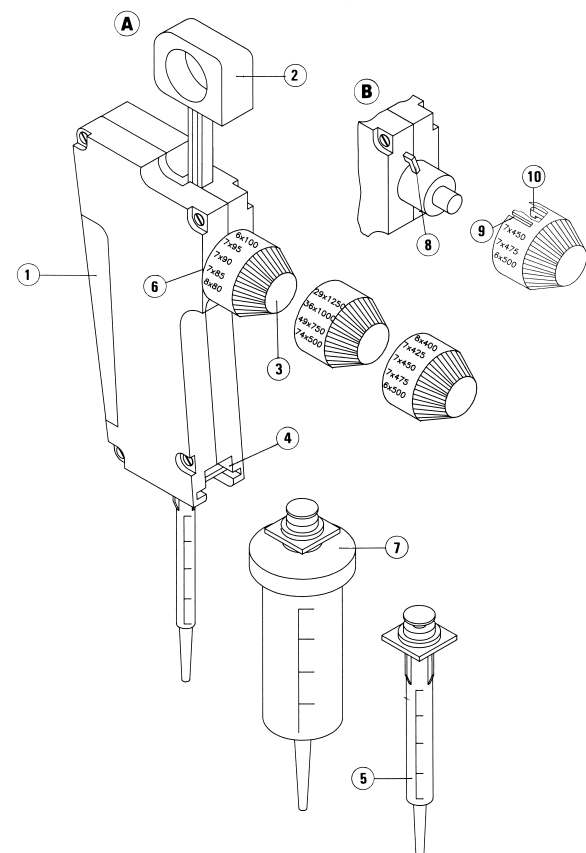
ACCESSOIRES

Description	Colour code	Volume	Cat. No.	Description	Packaging	Cat. No.
Replacement setting nob	Yellow	10 – 100 µL	411.582	Workstation	1/pk	320.411
	Blue	50 – 500 µL	411.583	Adaptor for red syringe	5/pk	1.411.7
	Red	500 – 5000 µL	411.584			

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